Genotyping Protocol: MMRRC 41409

Assay Type: PCR; can distinguish homozygous animals from heterozygous samples.

DNA Extraction: DNA from tail snips was extracted using the following protocol:

- 1. Place a ~3mm piece of mouse or rat tail in 180ul of 50mM NaOH in a 0.2mL tube and vortex
- 2. Incubate at 95°C for 10 minutes, then at 40°C for 5 minutes to prevent popped tube caps.
- Add 20ul of 1M Tris-HCl (pH 8.0) and vortex well. Use this lysate as your DNA template in KOD Xtreme PCR reactions.

Strain Description: This strain was created using a gene trap protocol. The vector pGT0lxf was inserted into intron 13 of mouse *Svil*. This vector contains an intron from the murine engrailed 2 gene upstream of a Beta-galactosidase/neomycin resistance (Beta-geo) fusion protein. Details can be found in Edelstein et al (2012) Circulation 125:2762-2771.

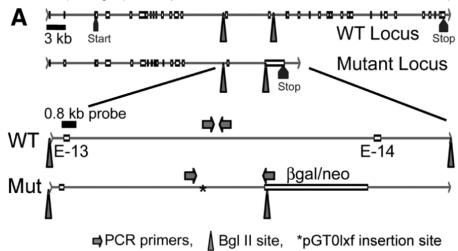


Image from Edelstein et al (2012) Circulation 125:2762-2771.

Primer Information:

1) Name: M41409 WT F
2) Name: M41409 WT R
3) Name: M41409 Mut F
4) Name: M41409 Mut R
Sequence: 5'-CAT GCT GAT GGA TTA GTG TTC TCA GGG ATA AGT CCT TTA C-3'
Sequence: 5'-GCA GAA GCA GAA TAC TGG CAG AAG GTC TAC TTT ACA TCT T-3'
Sequence: 5'-GCC TTT GGT ATT CTA CCC AAA AGA TGG TTA TAC CCA G-3'
Sequence: 5'-CAT CAA CAT TAA ATG TGA GCG AGT AAC AAC CCG TC-3'

Primer location: M41409 WT F and R and M41409 Mut F are located in Intron 14 of mouse *Svil.* M41409 Mut R is located in the inserted Beta galactosidase.

Assay Name: Svil KO PCR

Wild Type PCR Master Mix Components:

| component | manufacturer | concentration | μ l/rxn |
|-------------------|--------------------------|---------------|----------------|
| KOD Xtreme Buffer | Millipore | 2X | 10 |
| KOD Xtreme dNTPs | Millipore | 2mM | 4 |
| M41409 WT F | Sigma or IDT | 25µM | 0.3 |
| M41409 WT R | Sigma or IDT | 25µM | 0.3 |
| KOD Xtreme Taq | Millipore (Cat# 71975-3) | 1 U/μl | 0.4 |
| sterile water | | | 3 |

Mutant PCR Master Mix Components:

| component | manufacturer | concentration | μ l/rxn |
|-------------------|--------------------------|---------------|----------------|
| KOD Xtreme Buffer | Millipore | 2X | 10 |
| KOD Xtreme dNTPs | Millipore | 2mM | 4 |
| M41409 Mut F | Sigma or IDT | 25µM | 0.3 |
| M41409 Mut R | Sigma or IDT | 25µM | 0.3 |
| KOD Xtreme Taq | Millipore (Cat# 71975-3) | 1 U/μl | 0.4 |
| sterile water | | | 3 |

PCR Setup:

Final Reaction: 18 µl master mix & 2 µl extracted DNA (10-20ng DNA).

All reactions were performed in 200 μ l thin walled PCR tubes and were run in Perkin Elmer 2400 thermocycler or Applied Biosystems 2700 thermocycler.

Wild Type PCR Cycle Parameters:

| 1) | 95°C | 4 minutes |
|----|------|-----------|
| 2) | 94°C | 1 minute |
| 3) | 70°C | 1 minute |

4) 72°C 1 minute, 30 seconds
 5) Repeat steps 2-4 34 times for a total of 35 cycles

6) 72°C 7 minutes

7) 4°C hold until refrigerate product

Mutant PCR Cycle Parameters:

1) 95°C 4 minutes 2) 94°C 1 minute 3) 65°C 1 minute

4) 72°C 2 minutes, 30 seconds 5) Repeat steps 2-4 34 times for a total of 35 cycles

6) 72°C 7 minutes

7) 4°C hold until refrigerate product

Product Analysis:

All products were analyzed on the Qiaxcel (instrument and all supplies from Qiagen) with the Qiaxcel DNA Screening Kit (Cat# 929004).

Alignment and Size Markers:

Wild Type PCR: QX Alignment Marker 15bp/3Kb (Cat# 929522) Size Marker: QX DNA Size Marker 100-3Kb (Cat# 929553)


Mutant PCR: QX Alignment Marker 15bp/5Kb (Cat# 929524) Size Marker: QX DNA Size Marker 250-4Kb (Cat# 929562)

Both PCRs use: Method: AM320 Injection: 10s at 5KV, Separation: 320s at 6KV


Expected products:

Wild Type PCR: 800bp product Mutant PCR: 2.5kb product

Example Gels:

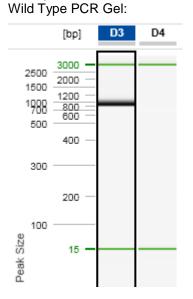

Mutant PCR Gel:

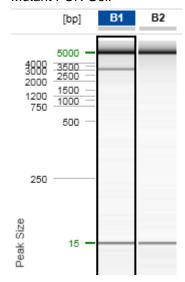

Lane D3 displays a sample positive for the Wild Type allele (800bp product)
Lane D4 displays a sample negative for the Wild Type allele (no product)

Please note: the 15bp and 3kb bands are reference markers specific to the QIAxcel method and do not represent expected products. Lane B1 displays a sample positive for the Mutant allele (2.5kb product)
Lane B2 displays a sample negative for the Mutant allele (no product)

Please note: the 15bp and 3kb bands are reference markers specific to the QIAxcel method and do not represent expected products.

Genotyping Analysis:

| Wild-type |
|-------------------|
| Heterozygous |
| Homozygous mutant |

Wild Type PCR 800bp product 800bp product no product

Mutant PCR no product 2.5kb product 2.5bp product